



In order to look up your data in the FACSuite software again later close the worklist and click in the worklist overview **Export > Worklist with data**.

Assay transfer

In case you would like to switch between FACSVerse and FACSLyric you can transfer your assays to the respective instrument. Choose in the Library  **Assays > User-Defined Assays**, click on the assay you want to transfer, then choose **File > Export** and save it in your folder on an external hard drive. On the other instrument choose in the Library **Assays > User-Defined Assays**, click onto the assay in your folder on the external hard drive, then choose **File > Import** and open it. Then you need to update your settings via the Setup&QC tab . Choose **Assay/Tube settings Setup**, select the correct bead lot and your assay, put a tube with FACSuite CS&T beads on the manual port and click on Run. Afterwards you can use your assay.

Questions and Problems

- ✓ Please call the FlowKon staff -3947/-3949/-2187 or write an email to flowkon@uni-konstanz.de when something is missing or not working properly.
- ✓ If an incident occurs during a session, please report it immediately in the PPMS incident sections so every user is aware of this and can re-plan his/her experiments if necessary.